



Midisart® 2000

The ready-to-use filter
for sterilizing gases and
venting

Midisart® 2000.

The sterilizing filter with impressive performance and total handling ease.



Midisart 2000 filtration units are ideal tools in biotechnology, the pharmaceutical industry, research institutes and anywhere you need sterile vents, bioisolation or sterile air and gases.

Midisarts are excellent for

- sterile venting of filling vessels and fermentation carboys, including culture vessels and CO₂ incubators (6 to 120 liters)
- venting of holding tanks for sterile, distilled water and liquid culture media
- autoclave venting
- in-line sterilization of and particulate removal from air and gases, such as sterilization of air for small fermenters

Midisart 2000 filtration units have been specially designed for maximum handling ease and safety. Tapered hose barbs ensure a simple and secure hold for 6- to 12-mm inner diameter tubing. Midisart is lightweight – only 20 g – so it will not weigh down or kink tubing.

User benefits:

1. Maximum handling ease

- Midisart 2000 comes individually packaged and presterilized – it's ready to connect!

2. Extra reliability and safety

- Midisart 2000 is integrity testable and delivers reproducible results.
- The membrane is reinforced with polypropylene gauze, giving the Midisart unit added stability and making it pressure resistant up to 3 bar (approx. 44 psi).
- Midisart 2000 entirely eliminates moisture breakthrough because of its inherently hydrophobic PTFE material.
- In addition, Midisart is biosafe because all materials of construction meet the requirements of the current USP Plastics Class VI testing.
- Midisart 2000 units easily withstand at least 20 autoclaving cycles with no loss in performance. The convenient Memory Discs supplied with each Midisart 2000 in UPN-coded boxes enable you to keep track of the number of autoclaving cycles by marking or clipping off each cycle. This feature is key in complying with GLP and ISO standards for traceable documentation.

3. Quality control certificate

- Each unit is automatically tested 100% for housing and membrane sealing during manufacture as part of our zero-defect quality control testing.
- The lot number and the individual unit number are imprinted on the top part of each Midisart 2000 housing to ensure complete traceability.

Midisart 2000 units are visually inspected before they are packaged.

In addition to 100% leak testing, random samples taken from each lot undergo the following tests to assure compliance with Sartorius' stringent in-house quality assurance standards:

- Housing burst pressure test
- Pressure-hold test
- Bubble point test
- Pyrogen test
- Sterile filtration capability
- Flow rate test
- Sterility test

Performance

- With a diameter measuring just 64 mm, Midisart incorporates a filter area of 20 cm², which means that it is "packed" with high flow rate performance power!
- Midisarts multiply filtration performance in more ways than one. They can be autoclaved at least 20 times at 134°C!

Chemical compatibility

The materials used in Midisart (PTFE and polypropylene) give it excellent compatibility with the solvents and other chemicals listed on the right. However, its compatibility can be affected by various factors, such as temperature, concentration, composition, etc. We therefore recommend that you perform a trial filtration run to test whether Midisart is compatible with the particular medium you wish to filter.

Midisart 2000 can also be used to filter aqueous solutions. In this case, it must be first wetted with alcohol to overcome the membrane's hydrophobicity.

- Acetic acid (concentrated), acetone, acetonitrile
- n-butanol
- Cellosolve (ethyl), chloroform
- Diethylacetamide, dimethyl formamide, dimethyl sulfoxide, dioxane
- Ethanol, ethyl acetate, ethylene glycol
- Freon TF
- Gasoline
- 1 N hydrochloric acid, hexane
- Isobutanol, isopropanol
- Methanol, methylene chloride, methyl ethyl ether, methyl ethyl ketone
- Sodium hydroxide (5%)
- Pentane
- Tetrahydrofuran, toluene, trichloroacetic acid, trichloroethane
- Water
- Xylene



Technical specifications

Filter material	PTFE – reinforced with polypropylene gauze	
Housing material	Polypropylene	
Filtration area	20 cm ²	
Housing diameter	64 mm	
Priming volume	Approx. 3 ml	
Maximum operating pressure	300 kPa (3 bar = 44 psi)	
Water penetration point (breakthrough)	0.2 µm – approx. 400 kPa (4 bar = 58 psi) 0.45 µm – approx. 300 kPa (3 bar = ~ 44 psi)	
Max. autoclaving temperature	134°C	
Max. autoclave cycles	20	
Hold-up volume	Before the bubble point	approx. 1.0 ml
	After the bubble point	approx. 0.5 ml
Biosafety	USP Plastics Class VI	
Bubble point with isopropanol (60%)	0.45 µm	≥ 0.9 bar (~13.1 psi)
	0.2 µm	≥ 1.1 bar (~16 psi)
Flow rate for air at Δp = 0.1 bar (1.45 psi) (1 bar = 100 kPa = 14.5 psi)	0.2 µm pore size	5.0 l/min
	0.45 µm pore size	8.5 l/min

Order information

Order numbers	Pore size	Membrane	Connectors E A	Pieces/Case	Sterile
17804 E	0.45 µm	PTFE	Olive Olive	12	Yes
17804 G	0.45 µm	PTFE	Olive Olive	25	Yes
17804 NPE	0.45 µm	PTFE	1/8" 1/8" NPT	12	Yes
17804 NPG	0.45 µm	PTFE	1/8" 1/8" NPT	25	Yes
17839 UNN	0.45 µm	PTFE	Luer Lock female Olive	100	No
17805 E	0.2 µm	PTFE	Olive Olive	12	Yes
17805 G	0.2 µm	PTFE	Olive Olive	25	Yes
17805 NPE	0.2 µm	PTFE	1/8" 1/8" NPT	12	Yes
17805 NPG	0.2 µm	PTFE	1/8" 1/8" NPT	25	Yes
17805 UPN	0.2 µm	PTFE	Olive Olive	100	No
17809 UNN	0.2 µm	PTFE	1/8" 1/8" NPT	100	No
17812 UNN	0.2 µm	PTFE	1/8" Olive	100	No

In the interest of further development of Sartorius products, we reserve the right to make changes without notice.

Sartorius AG
Weender Landstrasse 94–108
37075 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289

www.sartorius.com

Sartorius BBI Systems GmbH
Schwarzenberger Weg 73–79
34212 Melsungen, Germany

Phone +49.5661.71.3400
Fax +49.5661.71.3702

www.sartorius-bbi-systems.com

Sartorius North America Inc.
131 Heartland Blvd.
Edgewood, New York 11717, USA

Phone +1.631.254.4249
Toll-free +1.800.3687178
Fax +1.631.254.4253

Sartorius BBI Systems Inc.
999 Postal Road
Allentown, PA 18109, USA

Phone +1.610.2666262
Toll-free +1.800.258.9000
Fax +1.610.266.9319

Sartorius Ltd.
Longmead Business Park
Blenheim Road, Epsom
Surrey, KT19 9 QQ, U.K.

Phone +44.1372.737159
Fax +44.1372.726171

Sartorius S.A.
4, rue Emile Baudot
91127 Palaiseau Cedex, France

Phone +33.1.6919.2100
Fax +33.1.6920.0922

Sartorius S.p.A.
Via dell'Antella, 76/A
50011 Antella (FI), Italy

Phone +39.055.63.40.41
Fax +39.055.63.40.526

Sartorius, S.A.
C/Isabel Colbrand 10–12
Planta 4, Oficina 121
Polígono Industrial de Fuencarral
28050 Madrid, Spain

Phone +34.91.3586091
Fax +34.91.3588804

Sartorius Technologies N.V.
Luchthavenlaan 1–3
1800 Vilvoorde, Belgium

Phone +32.2.756.0670
Fax +32.2.756.0681

Sartorius K.K.
KY Building, 8–17
Kitashinagawa 1-chome
Shinagawa-ku
Tokyo 140-0001, Japan

Phone +81.3.3740.5407
Fax +81.3.3740.5406